

Olink®

COVID-19 1% TritonX-100 Inactivation protocol

Guidelines

Olink Proteomics, Analysis Service Uppsala and Analysis Service Boston accept COVID-19 patient samples for testing according to the following guidelines:

All samples should be TritonX-100 inactivated according to the following protocol PRIOR TO SHIPPING:

- Volume/sample for 1% TritonX-100 inactivation: 45 µL
- TritonX-100 and PBS
- Duration: 2 hours
- Temperature: Room temperature


NOTE: ALL samples that are being analyzed for a COVID-19 project should go through the same inactivation protocol, REGARDLESS of COVID-19 status (positive or negative).

Preparation

Overview

There are two steps in the protocol before shipping the COVID-19 samples to Olink:

1. Plating the samples according to Olink guidelines.
2. 1% TritonX-100 inactivation according to Olink guidelines.

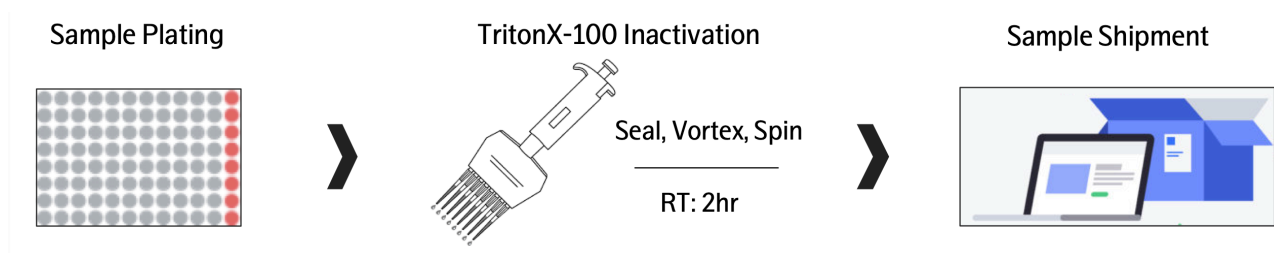


Required material and equipment

- 96-well PCR plate (compatible for up to 200 μL)
NOTE: Make sure the plate is -80°C , dry-ice resistant and easily resealable.
- Stock solution: 1 mL of 10% TritonX-100 in PBS per 96-well plate
- 8-well PCR strip
- Adhesive plastic film (temperature-resistant) **compatible with -80°C** , and heat resistant up to minimum $+56^{\circ}\text{C}$ (e.g. seal from Life Technologies #4306311)

Sample plating

- Pre-processing of blood samples: **Serum and plasma** need to be prepared according to standard protocols, aliquoted and inactivated before shipping to Olink. Whole blood or unprocessed serum/plasma samples cannot be sent.
- Plate **45 μL** of each sample in a temperature-resistant, non-protein binding plastics.
NOTE: Please note that the maximum sample volume recommended for shipment in plates is 100 μL .
- Ensure that your samples are randomized and balanced across all your plates in the order ready to be run and compatible with our plate layout (see below).
NOTE: For any assistance with randomization, please contact support@olink.com.

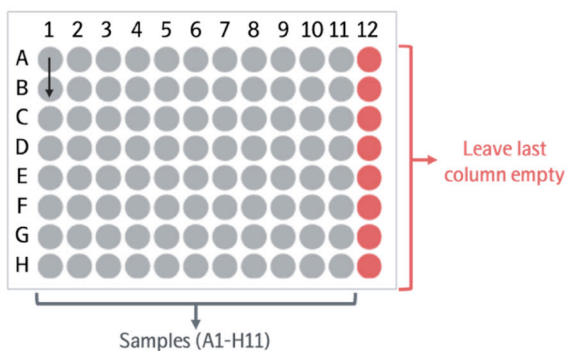

Protocol: 1% TritonX-100 inactivation

1. Prepare 10% TritonX-100 stock: Add 100 μL TritonX-100 into 900 μL PBS and pipette up and down slowly to mix.
2. Distribute 100 μL 10% TritonX-100 stock into each well of 8-well PCR strip (per sample plate).
3. To each sample well, add 10% TritonX-100 to final concentration of 1% using an 8-channel multichannel pipette:

Sample volume	10% TritonX-100 volume	Final volume
45 μL	5 μL	50 μL

4. Ensure that each well is separately sealed using an adhesive film or individual seals.
5. Vortex the plates thoroughly and spin down at 400g x 1min at RT.
6. Incubate the sample plate in the biosafety hood at RT for 2 hours.
7. Store your sample plates in -80°C freezer until ready for shipping.

Sample shipping

IMPORTANT: BEFORE SHIPPING THE SAMPLES, the COVID-19 protocol certification document needs to be signed and sent to Olink team. Please contact support@olink.com to receive this document.

- Samples should be shipped on dry ice, sufficient for shipment to the Analysis Service lab in Uppsala or Boston.
- Make sure the plates are securely sealed, add extra adsorption paper in case of leakage and make sure to place the sample plates in a sealed bag.
- Ensure that the heavy dry ice will not damage the samples during shipment.
- Contact your preferred courier for local regulations regarding appropriate labeling for risk assessment and shipment of your human samples on dry ice.
- Place a copy of the Work Order as well as the COVID-19 inactivation protocol certification in the parcel to identify samples.
- Send the tracking number and courier company name to the corresponding service unit: service@olink.com (Uppsala) or service_boston@olink.com (Boston), and add quote number in the communication.
NOTE: *Our business hours are Mon-Fri 8 a.m. to 5 p.m.*
- Upon arrival, the samples and documents will be examined, and you will receive a confirmation email.

www.olink.com

For research use only. Not for use in diagnostic procedures.

This product includes a license for non-commercial use. Commercial users may require additional licenses. Please contact Olink Proteomics AB for details.

There are no warranties, expressed or implied, which extend beyond this description. Olink Proteomics AB is not liable for property damage, personal injury, or economic loss caused by this product.

Olink® is a registered trademark of Olink Proteomics AB.

© 2020 Olink Proteomics AB. All third party trademarks are the property of their respective owners.

Olink Proteomics, Dag Hammarskjölds väg 52B, SE-752 37 Uppsala, Sweden

2020-09-17